

Create Account at traceablelive.com

1

Username


Password

[Forgot password?](#)

New Account Login

Fill out required fields

2

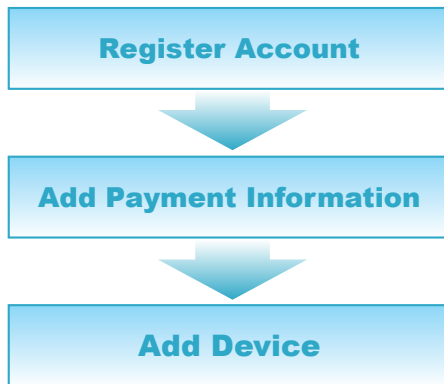

Back of unit

EXAMPLE 6500CC
S/N 100080013
KEY 5555123456
traceable.com 281-482-1714

By: 100 Due: Nov, 19 202

EXAMPLE

WiFi Setup

3




Press/HOLD WiFi button "AP" will appear on display.

Press WiFi button again, "AP UAT" will appear on display.

When "AP READY" appears, proceed to next step.

Use Internet enabled device to connect to WIFI

4

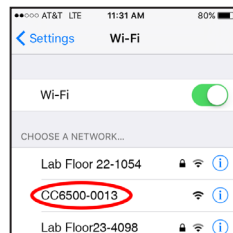

Select Network:
CC6500-0013
see example network

Example
Network: **CC6500-0013**

EXAMPLE 6500CC
S/N 100080013
KEY 5555123456
traceable.com 281-482-1714

By: 100 Due: Nov, 19 202


EXAMPLE

Open Browser, type
192.168.1.1 in address bar

5

Select local network and enter password.

Security type WPA2 recommended or
contact your network admin.

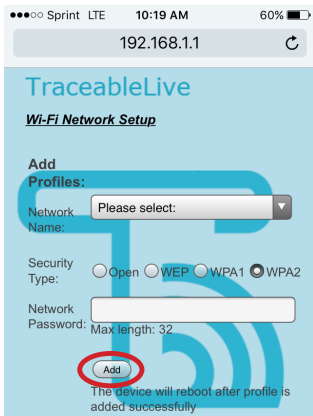

CONGRATULATIONS!
Your device will restart.

Still having trouble?
Please give us a call
281-482-1714 or email
support@traceable.com

TraceableLIVE®
QUICK START
GUIDE